



Presto FF-C18

250 x 4.6 mm


Application

**Mucopolysaccharides**


ムコ多糖

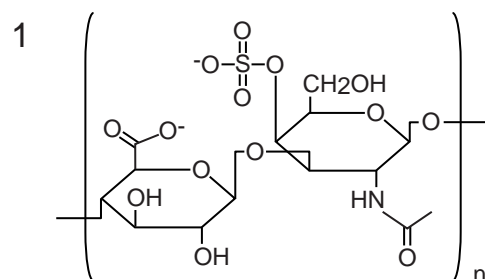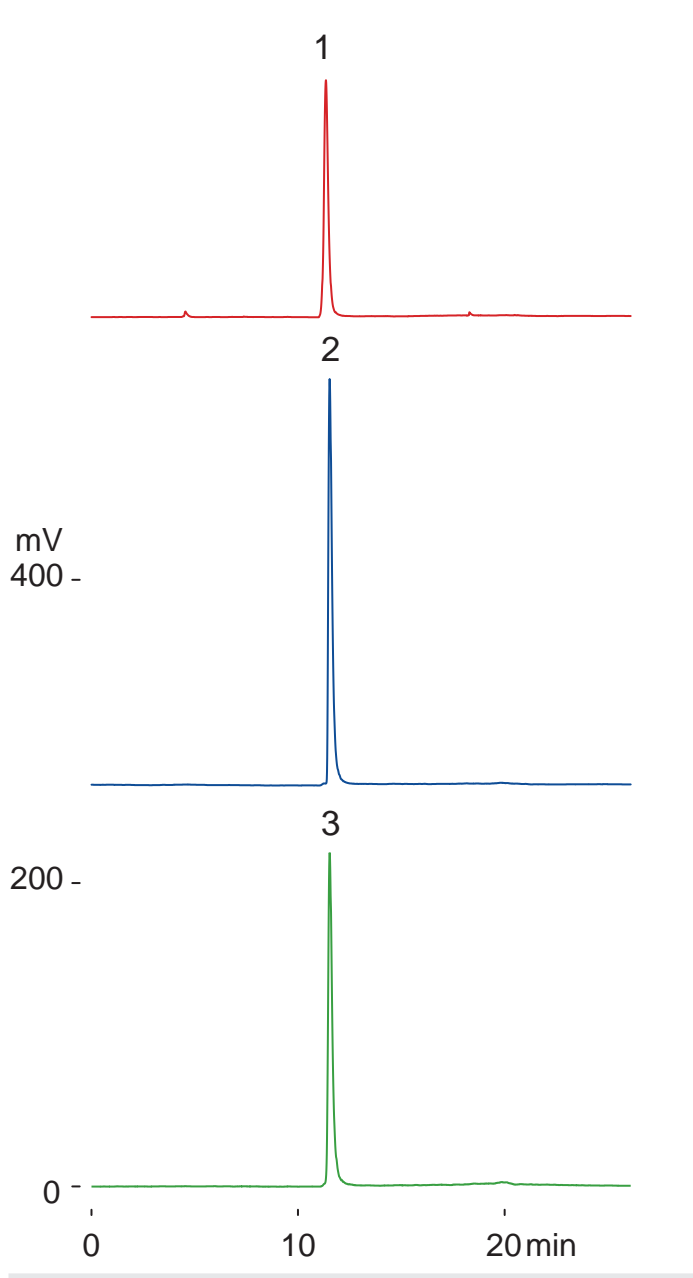

chondroitin sulfuric acid  
(~ 700 kDa)

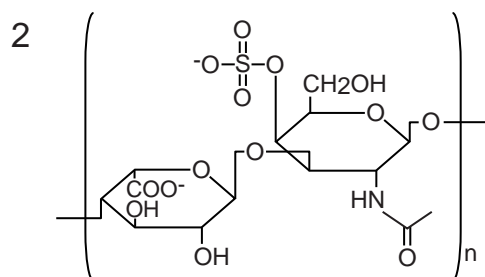

dermatan sulfuric acid  
(ca. 20 kDa)

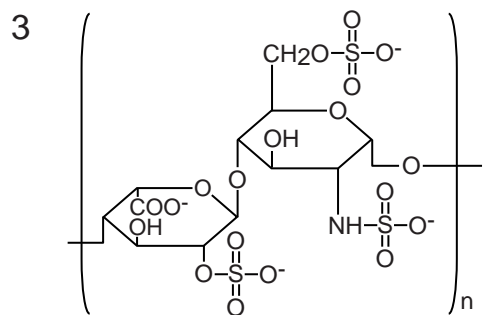

heparin (5 ~ 20 kDa)

Presto FF-C18, 250 x 4.6 mm

A: water /triethylamine /acetic acid = 100 /1.1 /0.5

B: acetonitrile

0-40 %B (0-15min)

0.35 mL/min (24 MPa), 37 deg.C

ELSD (spray chamber 50 deg.C, drift tube 100 deg.C)

2 uL (10 ug)