








Unison UK-Phenyl

100 x 3 mm


HPLC COLUMN APPLICATION

LC-MS: Neonicotinoid pesticides

ネオニコチノイド系殺虫剤の LC-MS 分析

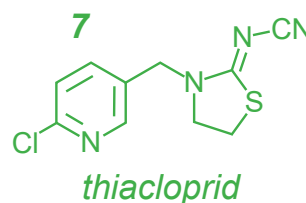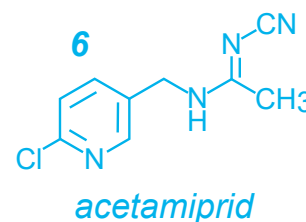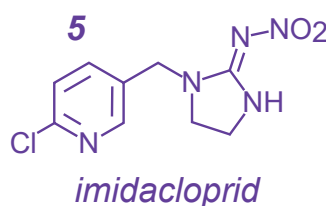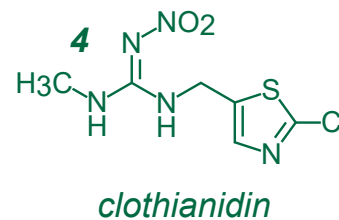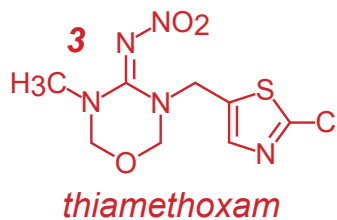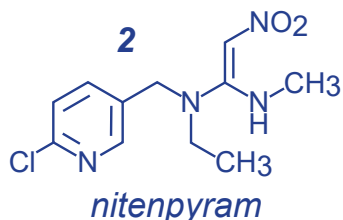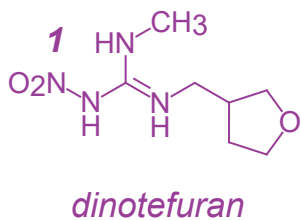

MZ-Analysentechnik GmbH
Barcelona-Allee 17 • D-55129 Mainz
Tel +49 6131 880 96-0
Fax +49 6131 880 96-20
e-mail: info@mz-at.de
www.mz-at.de

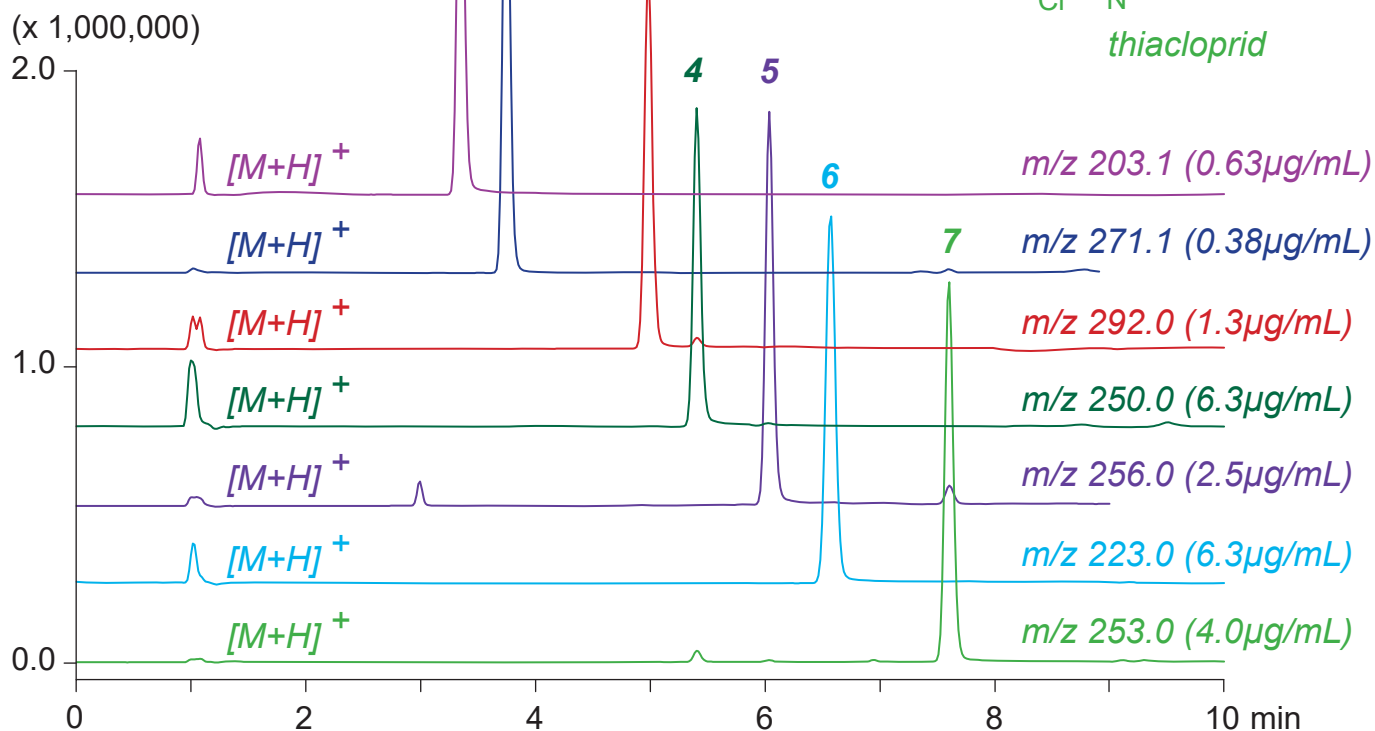

Unison UK-Phenyl, 100 x 3 mm
 A: water / formic acid = 100 / 0.1
 B: acetonitrile / formic acid = 100 / 0.1
 5-50%B (0-10min)
 0.4mL/min (6MPa), 37°C, Single Quad. MS (ESI)
 5µL (0.38-6.3µg/mL, 0.1% formic acid)