
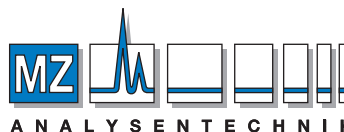


## UHPLC Analysis of Fat Soluble Vitamins on Titan™ C18, Acetonitrile Mobile Phase

column . . . . . TITAN C18, 50 mm x 2.1mm I.D. x 1.9 µm particles (577122-U)  
mobile phase . . . . . 100% acetonitrile  
flow rate . . . . . 0.4 mL/min  
pressure . . . . . 1755 psi (121 bar)  
column temp. . . . . 30 °C  
detector . . . . .  
injection . . . . . 0.5 µL  
sample . . . . . All compounds at 0.5 mg/mL except  
retinyl acetate at 0.05 mg/mL in methanol  
Application No. . . . . **G006221**

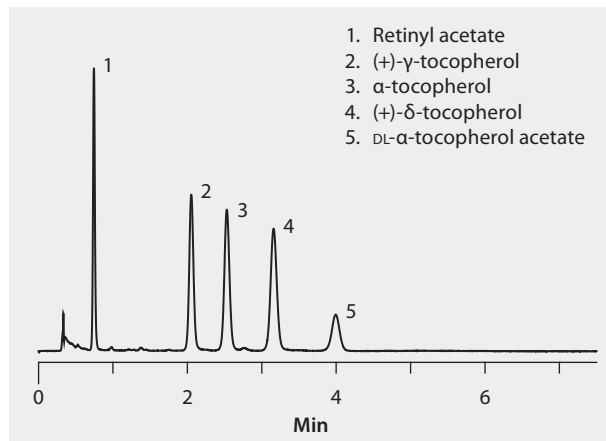

### AUTHORIZED DISTRIBUTOR

MZ-Analysentechnik GmbH, Barcelona-Allee 17 • D-55129 Mainz

Tel +49 6131 880 96-0, Fax +49 6131 880 96-20

e-mail: [info@mz-at.de](mailto:info@mz-at.de), [www.mz-at.de](http://www.mz-at.de)