



# EPA 8327 PFAS on Raptor C18

• Short cycle time (8 minutes).

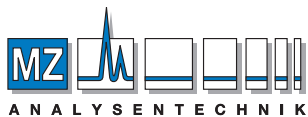

AUTHORIZED DISTRIBUTOR

MZ-Analysentechnik GmbH  
 Barcelona-Allee 17 • D-55129 Mainz  
 Tel +49 6131 880 96-0  
 Fax +49 6131 880 96-20  
 e-mail: info@mz-at.de  
 www.mz-at.de

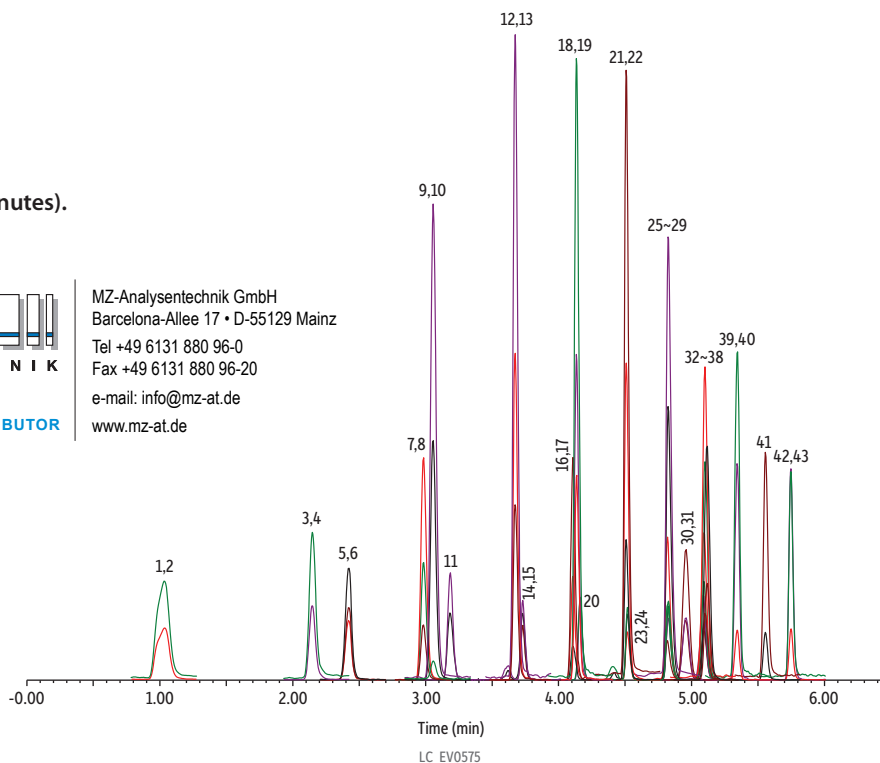

Peaks	tr (min)	Conc. (ng/L)	Precursor Ion	Product Ion	Product Ion	Column	Diluent	Conc.:	Inj. Vol.:	Mobile Phase	Time (min)	Flow (mL/min)	%A	%B
1. Perfluorobutanoic acid (PFBA)	1.04	80	231.1	169.1	-	Raptor C18 (cat.# 9304A52)	50:50 Water:methanol, 0.1% acetic acid	80 ppt (40 ppt for surrogate or isotopically-labelled PFAS)	10 µL	A: Water, 5 mM ammonium acetate B: Methanol	0.00	0.4	80	20
2. Perfluoro- <i>n</i> -[ <sup>13</sup> C <sub>4</sub> ]butanoic acid (M4PFBA)	1.04	40	217.0	172.0	-						6.00	0.4	5	95
3. Perfluoropentanoic acid (PFPeA)	2.14	80	263.0	219.0	-						6.01	0.4	80	20
4. Perfluoro- <i>n</i> -[ <sup>13</sup> C <sub>5</sub> ]pentanoic acid (M5PFPeA)	2.14	40	268.0	223.0	-						8.00	0.4	80	20
5. Perfluorobutane sulfonic acid (PFBS)	2.42	80	299.1	80.0	99.0									
6. Perfluoro-1-[2,3,4- <sup>13</sup> C <sub>3</sub> ]butyl sulfonic acid (M3PFBS)	2.42	40	302.0	80.0	-									
7. 1H, 1H, 2H, 2H-perfluorohexane sulfonic acid (4:2 FTS)	2.98	80	327.1	307.1	80.8									
8. 1H, 1H, 2H, 2H-perfluoro-1-[1,2- <sup>13</sup> C <sub>2</sub> ]hexyl sulfonic acid (M2-4:2 FTS)	2.98	40	329.0	309.0	-									
9. Perfluorohexanoic acid (PFHxA)	3.06	80	313.1	269.1	119.1									
10. Perfluoro- <i>n</i> -[1,2,3,4,6- <sup>13</sup> C <sub>5</sub> ]hexanoic acid (M5PFHxA)	3.06	40	318.0	273.0	-									
11. Perfluoropentane sulfonic acid (PFPeS)	3.19	80	349.1	80.0	99.0									
12. Perfluoroheptanoic acid (PFHpA)	3.67	80	363.2	319.1	169.1									
13. Perfluoro- <i>n</i> -[1,2,3,4- <sup>13</sup> C <sub>4</sub> ]heptanoic acid (M4PFHpA)	3.67	40	367.0	322.0	-									
14. Perfluorohexane sulfonic acid (PFHxS)	3.73	80	399.1	80.0	99.0									
15. Perfluoro-1-[1,2,3- <sup>13</sup> C <sub>3</sub> ]hexyl sulfonic acid (M3PFHxS)	3.73	40	402.0	80.0	-									
16. 1H, 1H, 2H, 2H-perfluorooctane sulfonic acid (6:2 FTS)	4.11	80	427.2	407.2	80.7									
17. 1H, 1H, 2H, 2H-perfluoro-1-[1,2- <sup>13</sup> C <sub>2</sub> ]octyl sulfonic acid (M2-6:2 FTS)	4.11	40	429.0	409.0	-									
18. Perfluorooctanoic acid (PFOA)	4.13	80	412.9	369.0	169.0									
19. Perfluoro- <i>n</i> -[ <sup>13</sup> C <sub>8</sub> ]octanoic acid (M8PFOA)	4.13	40	421.0	376.0	-									
20. Perfluoroheptane sulfonic acid (PFHpS)	4.17	80	449.2	80.0	99.0									
21. Perfluorononanoic acid (PFNA)	4.51	80	463.2	419.2	219.1									
22. Perfluoro- <i>n</i> -[ <sup>13</sup> C <sub>9</sub> ]nonanoic acid (M9PFNA)	4.51	40	472.0	427.0	-									
23. Perfluorooctane sulfonic acid (PFOS)	4.52	80	499.2	80.0	99.0									
24. Perfluoro-1-[ <sup>13</sup> C <sub>8</sub> ]octyl sulfonic acid (M8PFOS)	4.52	40	506.8	80.0	-									
25. 1H, 1H, 2H, 2H-perfluorodecane sulfonic acid (8:2 FTS)	4.81	80	527.2	507.2	80.8									
26. 1H, 1H, 2H, 2H-perfluoro-1-[1,2- <sup>13</sup> C <sub>2</sub> ]decyl sulfonic acid (M2-8:2 FTS)	4.81	40	529.0	509.0	-									
27. Perfluorononane sulfonic acid (PFNS)	4.83	80	549.2	80.0	99.0									
28. Perfluorodecanoic acid (PFDA)	4.83	80	513.2	469.2	219.1									
29. Perfluoro- <i>n</i> -[1,2,3,4,5,6- <sup>13</sup> C <sub>6</sub> ]decanoic acid (M6PFDA)	4.83	40	518.9	473.9	-									
30. N-methylperfluoro-1-octanesulfonamidoacetic acid (N-MeFOSAA)	4.95	80	570.2	419.2	483.2									
31. N-methyl-d3-perfluoro-1-octanesulfonamidoacetic acid (d3-N-MeFOSAA)	4.95	40	572.9	418.9	-									
32. Perfluorodecane sulfonic acid (PFDS)	5.09	80	599.2	80.0	99.0									
33. N-ethylperfluoro-1-octanesulfonamidoacetic acid (N-EtFOSAA)	5.09	80	584.2	419.2	483.1									
34. N-ethyl-d5-perfluoro-1-catanesulfonamidoacetic acid (d5-N-EtFOSAA)	5.09	40	589.0	419.0	-									
35. Perfluoroundecanoic acid (PFUnA)	5.10	80	563.2	519.2	269.1									
36. Perfluoro- <i>n</i> -[1,2,3,4,5,6,7- <sup>13</sup> C <sub>7</sub> ]undecanoic acid (M7PFUnA)	5.10	40	569.9	524.9	-									
37. Perfluorooctane sulfonic acid (FOSA)	5.12	80	498.2	78.0	-									
38. Perfluoro-1-[ <sup>13</sup> C <sub>8</sub> ]octanesulfonamide (M8FOSA)	5.12	40	505.9	78.0	-									
39. Perfluorododecanoic acid (PFDoA)	5.34	80	613.2	569.2	169.1									
40. Perfluoro- <i>n</i> -[1,2- <sup>13</sup> C <sub>2</sub> ]dodecanoic acid (M2PFDoA)	5.34	40	614.8	569.9	-									
41. Perfluorotridecanoic acid (PFTriA)	5.55	80	663.2	619.2	169.1									
42. Perfluorotetradecanoic acid (PFTreA)	5.75	80	713.2	669.2	169.1									
43. Perfluoro- <i>n</i> -[1,2- <sup>13</sup> C <sub>2</sub> ]tetradecanoic acid (M2PFTreA)	5.75	40	714.8	669.7	-									



© 2021 Restek Corporation. All rights reserved.  
[www.restek.com](http://www.restek.com)

Restek patents and trademarks are the property of Restek Corporation. (See [www.restek.com/Patents-Trademarks](http://www.restek.com/Patents-Trademarks) for full list.) Other trademarks in Restek literature or on its website are the property of their respective owners. Restek registered trademarks are registered in the U.S. and may also be registered in other countries.