

# A Fast Efficient Method to Determine the Presence of Nitrosamines in Cosmetics, Personal Care, and Consumer Products

Michael Jones<sup>1</sup> and Christopher Glover<sup>2</sup>
<sup>1</sup>Waters Corporation, Manchester, UK
<sup>2</sup>Northumbria University, Newcastle upon Tyne, UK

#### **APPLICATION BENEFITS**

This application note describes a method for the identification and quantitation of carcinogenic nitrosamines in consumer products offering:

- Increased sample throughput and a reduction of solvent usage due to reduced run times.
- A single, direct LC-MS method to analyze both non-volatile and also volatile nitrosamines with no requirements for derivitization.
- Quantitation of N-nitrosamines below the regulatory limit of 50 μg/kg.
- The ability to mine for additional unknowns utilizing RADAR™ data acquisition.

# WATERS SOLUTIONS

ACQUITY UPLC® I-Class System

ACQUITY UPLC HSS T3 Column

Xevo® TO-S micro

MassLunx® MS Software

Oasis® HLB SPE Cartridges

#### **KEY WORDS**

Nitrosamines, personal care products, cosmetics, consumer products, regulated substances, RADAR, PICS

#### INTRODUCTION

N-nitrosamines are a class of compounds that have been shown to exhibit carcinogenic and mutagenic effects in animal models at several different tissue sites and by several different routes of exposure. Nitrosamines can potentially be formed in consumer products, cosmetics and personal care formulations, either during manufacture or product storage. N-Nitrosamine formation occurs when secondary amines are present in addition to a nitrating agent such as nitrous acid, nitrites, or nitrogen oxides, generally under acidic conditions. Analysis of cosmetic, personal care and consumer products for nitrosamine content should be included in Product Safety Assessments to ensure compliance with the new Cosmetic and Consumer Products Regulations. It has also been noted that cosmetics over five years old had higher concentrations of N-nitrosamines than new samples of the same products. This indicates that the formation of N-nitrosamines limits the shelf life of cosmetics and should be assessed as part of a cosmetic product's routine safety evaluation.

In both the United States and the European Union, consumer products manufacturers ensure product safety prior to commercialization, list all ingredients on the product label, and comply with any restrictions that are established for ingredients and products. Any potential risk from a product is assessed as part of its safety evaluation. In the European Union, EU Directive 1223/2009 has been in force since 2013.<sup>5</sup> It strengthens the safety of cosmetic products and streamlines the framework for all operators in the sector. It provides a robust, internationally recognized regime, which reinforces product safety while taking into consideration the latest technological developments. EU 1223/2009 specifies in Annexes II and III that N-nitrosamines (along with other contaminants) must not be present in cosmetic products above levels of  $50 \,\mu g/kg$ . In the United States, the U.S. FDA, under the auspices of The Food, Drug, and Cosmetics Act (1938) and The Fair Packaging and Labelling Act (1966) considers any cosmetic product containing N-nitrosamines as adulterated, and as such, has the power to apply for those products to be seized and removed from the market.<sup>7,8</sup> In Asia, the ASEAN nations have modeled the ASEAN Cosmetics Directive<sup>9</sup> on the above European legislation. China, India, and Japan also regulate cosmetics in a similar fashion.

1

# [APPLICATION NOTE]

EU Directive 2009/48/EC pertains to the safety of toys, finger paints, and elastomeric materials likely to be placed in the mouth of children. This legislation also limits the cumulative presence of N-nitrosamines to less than  $50 \, \mu g/kg$  in total. The legislation applies to any toys or equipment containing elastomeric materials such as soothers and bottle teats aimed at children. The applicability of this method was also tested against analysis of infant feeding bottle teats for the presence of N-nitrosamines utilising sample preparation procedures set out in draft European Standard prEN 71 - 12:2012:E. Previously utilised methods for the analysis of N-nitrosamines in both cosmetics formulations and elastomeric materials include GC-TEA, GC-MS, and HPLC coupled to colorimetric detectors and LC-ESI-MS. Texture times ranged from  $12 \, \text{to} 35 \, \text{mins}$  for LC methods and  $20 \, \text{to} 40 \, \text{mins}$  for GC methods. TEA analysis was reported to be very time consuming, which limited sample throughput. It was also generally only applicable to analysis of volatile nitrosamines.

In this application note, we describe the analysis of eight volatile and non-volatile N-nitrosamines previously detected in cosmetics and personal care formulations using Waters® ACQUITY UPLC I-Class System coupled to Xevo TQ-S micro. The described method offers several benefits over previous methodologies including enhanced selectivity and sensitivity over colorimetric and ESI-MS methods. Quantitation is possible below the regulatory limits of  $50 \, \mu g/kg$ . Runtimes are reduced offering associated time and cost savings over HPLC and GC methods. The RADAR acquisition method used allows the ability to mine simultaneously acquired full scan data to search for other contaminants.

# **EXPERIMENTAL**

Cosmetic and personal care samples were prepared for analysis by an extraction protocol developed and validated by Qiang et al.<sup>11</sup> The samples underwent sonic extraction followed by sample cleanup utilizing Oasis HLB SPE Cartridges. Liquid and gel samples were prepared using 1:1 MEOH: acetone as a sonic extraction medium. Solid or semi-solid cosmetic samples were prepared using ethanol: acetone 4:1.

# Sonic extraction protocol

- 1g aliquot of a personal care formulation was dissolved in a 5.0 mL appropriate extraction medium with vortex mixing and sonicated for 30 min.
- The resulting extraction mixture was centrifuged at 6000 rcf for 15 min.
- 2-mL supernatant was evaporated to dryness under a constant stream of nitrogen at 40 °C, and reconstituted in 2 mL 30% aqueous methanol ready for solid phase extraction.

# SPE protocol

- A 6 cc<sup>3</sup> Oasis HLB SPE Cartridge (<u>Part no. WAT106202</u>)
   containing 200 mg sorbent was primed with 6 mL methanol,
   followed by 6 mL ultrapure water.
- The sample was loaded onto the SPE cartridge and washed with 3 mL 30% aqueous methanol. The sample was then extracted from the cartridge with 6 mL MeOH.
- The sample was prepared for injection by evaporating to dryness under N<sub>2</sub>, reconstituting in 1 mL 2.5% aqueous methanol and filtration through a 0.2 μm membrane filter.

# LC conditions

LC system:	ACQUITY UPLC I-Class
------------	----------------------

Run time: 9.1 min

Column: ACQUITY UPLC HSS T3

1.7 μm, 2.1 x 100 mm

Column temp.: 40 °C Sample temp.: 5 °C

Mobile phase A: Water (0.1% formic acid)

Mobile phase B: Methanol (0.1% formic acid)

20 µL

Flow rate: 0.3 mL/min

Injection volume:

Mobile phase gradient is detailed in Table 1.

	Time (min)	Flow rate (mL/min)	%A	%В	Curve
1	Initial	0.30	97.5	2.5	_
2	1.0	0.30	97.5	2.5	6
3	2.0	0.30	50.0	50.0	6
4	3.0	0.30	50.0	50.0	6
5	3.5	0.30	2.5	97.5	6
6	6.0	0.30	2.5	97.5	6
7	6.1	0.30	97.5	2.5	6

Table 1. ACQUITY I-Class mobile phase gradient.

#### MS conditions

Xevo TQ S-micro
APCI positive
15 V
150 °C
600 °C
1200 L/hr
200 L/hr
MRM with RADAR

Ionization parameters and transition pairs were optimized automatically using the IntelliStart™ function of MassLynx Software. IntelliStart is a feature within MassLynx that can be used to monitor system health, perform QC checks, and aid the method development process by increasing ease of use and decreasing user to user variation. In this instance the automatic tuning parameters feature of IntelliStart was used to determine optimum ionization parameters such as cone voltages for parent ions and collision gas energies for MRM transitions. Table 2 summarizes the optimum parameters for the analytes tested, along with information about the analyte names, chemical formulas, and CAS numbers.

Name, formula, CAS no.	Structure	Retention time (min)	Cone voltage (V)	Transition	Collision energy
N-Nitroso-dimethylamine, $C_2H_6N_2O$ , 62-75-9	N-O N	2.1	30	75.0→58.0*	13
N-Nitroso-diethylmine,	N O	3.1	24	103.1→47.0	15
C <sub>4</sub> H <sub>12</sub> N <sub>2</sub> O, 55-18-5		3.1	24	103.1→75.0*	9
N-Nitroso-dipropylamine,	N <sup>-,O</sup>	4.2	20	131.1→43.1	11
C <sub>6</sub> H <sub>14</sub> N <sub>2</sub> O, 621-64-7	/\/\\\	4.2	26	131.1→89.1*	8
N-Nitroso-diethanolamine,	Ν̈́O	1.4	10	135.0→74.0	8
$C_4 H_{10} N_2 O_3$ , 1116-54-7	но \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \	1.4	10	135.0→104.0*	4
N-Nitroso-dibutylamine,	ŃΣO	4.5	26	159.1→57.1*	12
C <sub>8</sub> H <sub>18</sub> N <sub>2</sub> O, 924-16-3	\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\\	4.5	20	159.1→103.0	9
N-Nitroso-pyrrolidine,	N = O	2.8	44	101.0→40.5	19
C <sub>4</sub> H <sub>8</sub> N <sub>2</sub> 0, 930-55-2	$\langle n \rangle$	2.0	44	101.0→55.0*	13
N-Nitroso-piperidine,	N.O	3.2	36	115.1→40.9	25
C <sub>5</sub> H <sub>10</sub> N <sub>2</sub> 0, 100-75-4		3.2	30	115.1→69.1*	17
N-Nitroso-morpholine,	N <sup>2O</sup>	2.6	28	117.0→45.0*	13
C <sub>4</sub> H <sub>8</sub> N <sub>2</sub> O <sub>2</sub> , 59-89-2		2.0	۷٥	117.0→85.9	9

Table 2. Aliphatic, cyclic, and heterocyclic N-nitrosamines previously found in consumer product formulations, structures, associated CAS numbers and empirical formulas, expected retention times, cone voltages, MRM transitions, and associated collision energy values.

# Instrument control, data acquisition, and results processing

MassLynx Software was used to control the ACQUITY UPLC I-Class System and the Xevo TQS-micro, and also for data acquisition. Data quantitation was achieved using TargetLynx™ Application Manager.

<sup>\*</sup>Refers to the quantitation transition.

#### RESULTS AND DISCUSSION

The analysis of eight N-nitrosamines historically found in cosmetic formulations was achieved using the Xevo TQ S-micro in MRM mode with APCI ionization in positive mode, coupled to an ACQUITY I-Class System.

The MRM transitions for each of the analytes were optimized using the IntelliStart function of MassLynx Software as previously discussed. LC conditions (column chemistry, gradient profile, injection volume, etc.) were based on previous work by Al-Kaseem et al<sup>16</sup>, and optimized to improve run time, elution order, and signal intensity.

Example chromatograms of standards prepared in the cosmetic matrix at a concentration equivalent to the regulatory limit of  $50 \,\mu g/kg$  are shown in Figure 1.

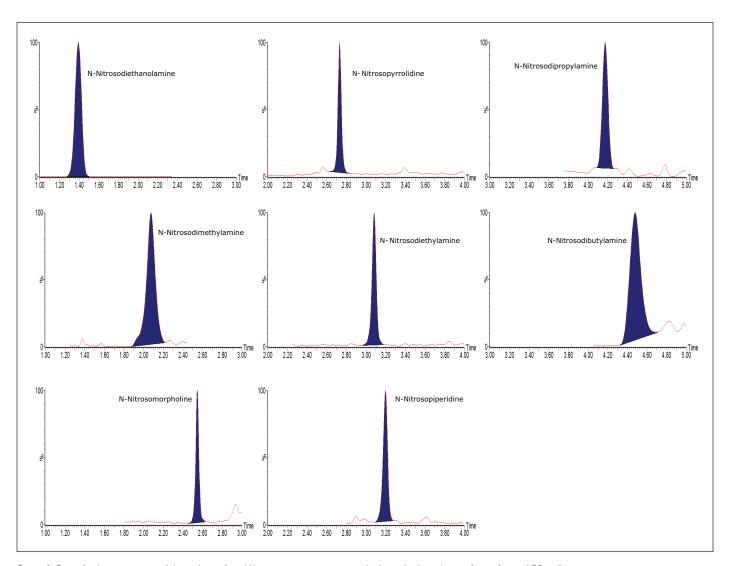


Figure 1. Example chromatograms of the eight analyzed N-nitrosamines in matrix spiked standards at the regulatory limit of 50  $\mu$ g/kg.

Mixed calibration standards from 0.001 ppm to 0.5 ppm were prepared and analyzed for all of the compounds considered. All nitrosamines were found to give a linear response over concentration ranges from 0.01 ppm to 0.5 ppm (10 times the regulatory limit of 0.05 ppm).  $R^2$  values for all curves were >0.99, and matrix effects were found to be <8%.

## Analysis of cosmetics and personal care samples

Samples were prepared from a range of cosmetic and personal care formulations including shampoo, hand wash, and lipstick. All samples were tested for the presence of the N-nitrosamines as detailed in the Experimental section, and were found to be free of N-nitrosamines. In order to test the suitability of the method, blank matrix samples were spiked with nitrosamines in amounts at or around the regulatory limit. In the example shown below in Figure 2, a shampoo sample was prepared, and spiked with N-nitrosodipropylamine at a concentration of 0.06 ppm (ng/mL). After sonic extraction and SPE the sample was analyzed alongside the matrix standards using a TargetLynx method to determine the N-nitrosodipropylamine concentration.

The data was quantified using the TargetLynx method to determine the amount of N-nitrosodipropylamine present in the spiked sample. Figure 1 shows the results generated in TargetLynx for the spiked shampoo sample, along with the calibration curve generated from the matrix standards to quantify the spiked sample. Figure 1 also shows the MRM trace for the confirmatory transition for N-nitrosodipropylamine. All of the results generated from the method for spiked samples were accurate in the range of 98% to 102% expected result for tested formulations including shampoo, hand wash, and lipstick.

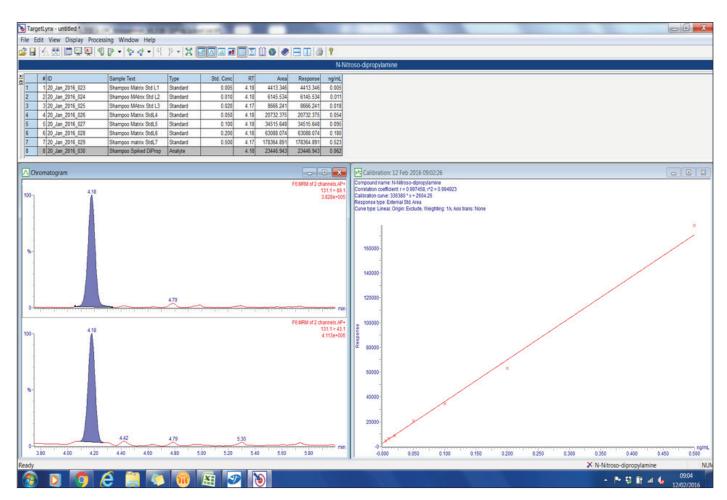


Figure 2. TargetLynx Result Summary for a shampoo sample spiked with N-nitrosodipropylamine.

## Analysis of baby feeding bottle teats

Silicon-free feeding teats from a leading manufacturer were tested according to a draft protocol set out in European Standard prEN 71 - 12:2012:E. 10 cm<sup>2</sup> portions of the elastomeric material used to form the teat were submerged in synthetic saliva containing a source of nitrite. Nitrite has been proven to convert secondary amines to N-nitrosamines under acidic conditions such as those found in the human stomach.<sup>17</sup> The teat samples were submerged for 30 minutes before the resultant supernatant was analyzed according to the method with no further sample preparation.

Unadulterated submerged samples were analyzed to determine the presence of nitrosamines. The samples were found to contain trace amounts of N-nitrosodibutylamine at levels approximately 10 times below the regulatory limits. In order to further test the efficacy of the method, additional teat samples were prepared and spiked with N-nitrosodiethanolamine at a concentration 1.4 times the regulatory limit. Samples were also prepared spiked with a mixture of three different N-nitrosamines, each below the regulatory limit, but with a cumulative concentration above the regulatory limit of  $50 \,\mu\text{g/kg}$ . Figure 3 shows The TargetLynx Result Summary for the N-nitrosodiethanolamine spiked sample, comprised of the calculated injections, a chromatogram of the spiked sample, and the N-nitrosodiethanolamine calibration curve generated for the calculation.

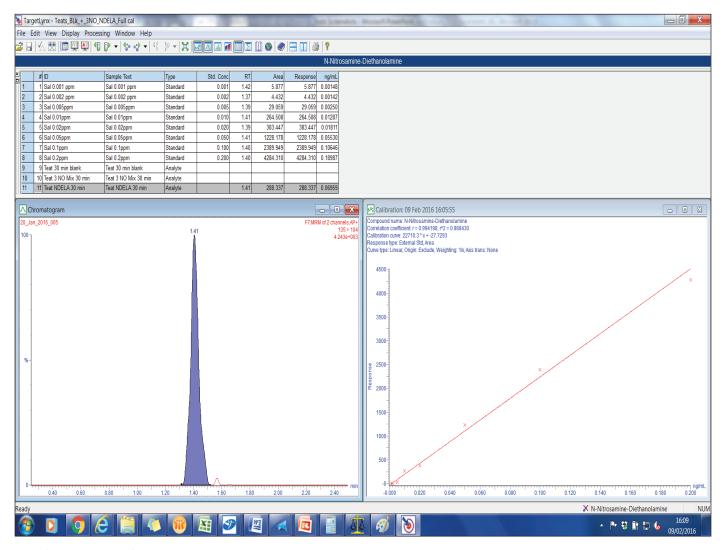


Figure 3. TargetLynx Result Summary for NDELA spiked baby bottle feeding teat sample.

The TargetLynx Sample Quantification reports generated from the NDELA spiked samples and the low level spiked samples are shown in Figure 4.

# Name	Trace	RT	Area	IS Area	Response	ng/mL	
N-Nitrosodimethyl amine	75 > 58						
2 N-Nitroso-diethylamine	103.1 > 75	3.09	576.317		576.317	0.018	
N-Nitroso-dipropylamine	131.1 > 89.1	4.19	1685.821		1685.821	0.022	
N-Nitroso-dibutylamine	159.1 > 57.1	4.54	3012.368		3012.368	0.008	
N-Nitrosamine-Diethanolami	. 135 > 104						
N-Nitroso-morpholine	117 > 45						
N-Nitroso-piperidine	115.1 > 69.1	3.21	938.489		938.489	0.017	
8 N-Nitroso-pyrollidine	101.1 > 55						
ama: 20 Jan 2016 005	Date: 20 Jan 2	016 Time:	40:34:44 ID	· Toot NDE	I A 30 min	Descripti	ion: Toot NDEL A 30 min
							ion: Teat NDELA 30 min
	Trace	016, Time:	10:34:44, ID	: Teat NDE	LA 30 min,	Descripti	ion: Teat NDELA 30 min
# Name							ion: Teat NDELA 30 min
# Name 1 N-Nitrosodimethyl amine	Trace						ion: Teat NDELA 30 min
# Name 1 N-Nitrosodimethyl amine 2 N-Nitroso-diethylamine	Trace 75 > 58						ion: Teat NDELA 30 min
# Name 1 N-Nitrosodimethyl amine 2 N-Nitroso-diethylamine 3 N-Nitroso-dipropylamine	Trace 75 > 58 103.1 > 75						ion: Teat NDELA 30 min
# Name  1 N-Nitrosodimethyl amine  2 N-Nitroso-diethylamine  3 N-Nitroso-dipropylamine  4 N-Nitroso-dibutylamine	Trace 75 > 58 103.1 > 75 131.1 > 89.1 159.1 > 57.1	RT	Area		Response	ng/mL	ion: Teat NDELA 30 min
# Name  1 N-Nitrosodimethyl amine  2 N-Nitroso-diethylamine  3 N-Nitroso-dipropylamine  4 N-Nitroso-dibutylamine  5 N-Nitrosamine-Diethanolami	Trace 75 > 58 103.1 > 75 131.1 > 89.1 159.1 > 57.1	RT 4.54	Area 2485.766		Response 2485.766	ng/mL 0.006	ion: Teat NDELA 30 min
# Name  1 N-Nitrosodimethyl amine  2 N-Nitroso-diethylamine  3 N-Nitroso-dipropylamine  4 N-Nitroso-dibutylamine  5 N-Nitrosomine-Diethanolami  6 N-Nitroso-morpholine  7 N-Nitroso-piperidine	Trace 75 > 58 103.1 > 75 131.1 > 89.1 159.1 > 57.1 135 > 104	RT 4.54	Area 2485.766		Response 2485.766	ng/mL 0.006	ion: Teat NDELA 30 min

Figure 4. TargetLynx Summary reports detailing the presence of N-nitrosamines in baby bottle feeding teats.

# RADAR analysis of aged consumer goods samples

As mentioned previously, it has been noted that N-nitrosamine levels can increase in personal care formulations upon storage.<sup>6</sup> The ability to determine the emergence of any N-nitrosamines over time would allow manufacturers to determine safe shelf lives for their products. To this end the method was used to test for the presence of nitrosamines in a new lipstick formulation and an old sample of the same brand. In addition to the targeted searching of the eight nitrosamines outlined in the method, utilizing the RADAR functionality available on the Xevo TQ-S micro instrument allows for simultaneous recording of multiple MRM transitions, along with scanning experiments across a defined *m/z* range. Figure 5 shows the TIC chromatograms comparing the new and aged lipstick samples. There appears to be an additional peak in the TIC with a retention time of approximately 4.55 minutes that does not correspond to N-nitrosodibutylamine, as shown by the absence of any peak at 4.55 minutes in the MRM transition chromatogram for that compound.

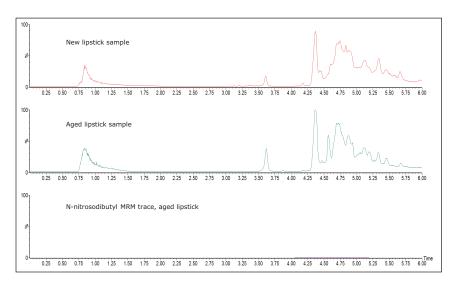


Figure 5. TIC chromatograms of new and aged lipstick formulation samples and an MRM transition chromatogram of N-nitrosodibutylamine showing its absence in the aged sample.

# [APPLICATION NOTE]

When analyses are conducted using RADAR data acquisition mode, expected analytes can be accurately quantified while simultaneously probing the sample for the presence of unknowns — without any loss of data integrity. Figure 6 again shows the TIC chromatograms for both the new and aged lipstick samples. When run in RADAR mode, the data acquired allows for examination of the relevant mass spectra to determine any differences at the observed peak time. The magnified MS spectra views in Figure 6 show the significant presence of an ion at m/z 199 in the aged sample over the new sample. Extracting this ion gives the extracted ion chromatogram shown in Figure 6 with a corresponding retention time of 4.55 minutes. A search of the literature suggests that m/z 199 possibly corresponds to the [M+H] $^+$  ion for N-nitrosodiphenylamine if the peak appearing in aged samples is due to N-nitrosamine formation. Further work is planned utilizing high resolution mass spectrometry (HRMS) to confirm the identity of the unknown ion present in the aged lipstick samples.

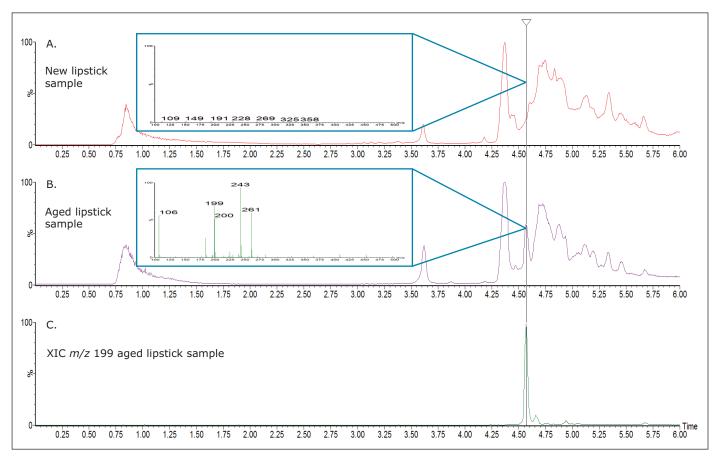


Figure 6. Probing the presence of unknowns using RADAR data acquisition: A. TIC of new lipstick sample and averaged MS spectra at 4.5 to 4.55 min.; B. TIC of new lipstick sample and averaged MS spectra at 4.5 to 4.55 min.; C. XIC of ion at m/z 199 in the aged lipstick sample.

# [APPLICATION NOTE]

#### CONCLUSIONS

An LC-APCI-MS method has been developed to rapidly detect and quantify eight commonly found nitrosamines in consumer product formulations. All analytes are quantifiable at or below the regulatory limits of  $50~\mu g/kg$  for cosmetics formulations and elastomeric toys. Three cosmetic formulations and one elastomeric material sample were tested for the presence of nitrosamines. The shampoo and hand wash formulations tested were found to be free of nitrosamines. The tested baby feeding teat sample was found to contain trace amounts of an N-nitrosamine below the regulatory limit. The incorporation of RADAR data acquisition into the method allows for the mining of full scan MS data for the potential presence of additional nitrosamines and suggested the presence of an additional N-Nitrosamine in the aged cosmetic formulation.

#### References

- 1. Ashley L M, Halver J E. Dimethylnitrosamine-induced hepatic cell carcinoma in rainbow trout. *J National Cancer Institute.* 41 (2): 531–552, 1968.
- 2. Barnes J M, Magee P N. Some Toxic Properties of Dimethylnitrosamine. British Journal of Industrial Medicine. 11 (3): 167–174, 1954.
- Jakszyn P, González C A. Nitrosamine and related food intake and gastric and oesophageal cancer risk: a systematic review of the epidemiological evidence. World Journal of Gastroenterology. 12 (27): 4296, 2006.
- Douglass M, Kabacoff B, Anderson G, Cheng M. The chemistry of nitrosamine formation, inhibition and destruction. J Soc Cosmet Chem. 29 (9): 581–606, 1978.
- Regulation (EC) No 1223/2009 of the European Parliament. Official Journal of the European Union, L342 (22.12.2009), 59–209, 2009.
- Matyska M T, Pesek J J, Yang L. Screening method for determining the presence of N-nitrosodiethanolamine in cosmetics by open-tubular capillary electrochromatography. *Journal of Chromatography A*. 887 (1): 497–503, 2000.
- 7. Administration, U. S. F. D., Fed Reg 44. 1979, 21365.
- 8. Administration, U. S. F. D., Fed Reg 58. 1983, 28288–28292.
- $9. \ \ \, \underline{\text{http://aseancosmetics.org/docdocs/directive.pdf}}$
- 10. Directive 2009/48/EC of the European Parliament and of the Council of 18 June 2009 on the safety of toys. Official Journal Journal of The European Union. L170 (30.6.2009): 1–37, 2009.
- Ma Q, Xi H W, Wang C, Bai H, Xi G C, Su N, Xu L Y, Wang J B. Determination of Ten Volatile Nitrosamines in Cosmetics by Gas Chromatography Tandem Mass Spectrometry. Chinese Journal of Analytical Chemistry. 39 (8): 1201–1207, 2011.

- 12. Flower C, Carter S, Earls A, Fowler R, Hewlins S, Lalljie S, Lefebvre M, Mavro J, Small D, Volpe N. A method for the determination of N-nitrosodiethanolamine in personal care products collaboratively evaluated by the CTPA Nitrosamines Working Group. *Intl J Cosmetic Sci.* 28 (1): 21–33, 2009.
- Schothorst R, Somers H. Determination of N-nitrosodiethanolamine in cosmetic products by LC-MS/MS. Analytical and bioanalytical chemistry. 381 (3): 681–685, 2005.
- 14. Schothorst R C, Stephany R W. Occurrence of N-nitrosodiethanolamine (NDELA) in cosmetics from the Dutch market. *Intl J Cosmetic Sci.* 23 (2): 109–114, 2001.
- Stepanov I, Carmella S G, Hecht S S, Duca G. Analysis of Tobacco-Specific Nitrosamines in Moldovan Cigarette Tobacco. *J Agricultural and Food Chemistry*. 50 (10): 2793–2797, 2002.
- 16. Al-Kaseem M, Al-Assaf Z, Karabeet F O. A Rapid, Validated RP-HPLC Method for the Determination of Seven Volatile N-nitrosamines in Meat. *Pharmacology & Pharmacy*. 5(3): 11, 2014.
- 17. Krul C A, Zeilmaker M J, Schothorst R C, Havenaar R. Intragastric formation and modulation of N-nitrosodimethylamine in a dynamic in vitro gastrointestinal model under human physiological conditions. Food and Chemical Toxicology. 42 (1): 51–63, 2004.



THE SCIENCE OF WHAT'S POSSIBLE.

Waters, ACQUITY UPLC, Xevo, Oasis, MassLynx, and The Science of What's Possible are registered trademarks of Waters Corporation.

RADAR, IntelliStart, and TargetLynx are trademarks of Waters Corporation. All other trademarks are the property of their respective owners.

Waters Corporation 34 Maple Street Milford, MA 01757 U.S.A. T: 1 508 478 2000 F: 1 508 872 1990 www.waters.com